



Here are your new **Genesis Ultra TLD-BP Dosimeters** for the next wear period.
Follow the instructions below to remove your old Blister Pack (from the reusable holder) and insert your new one.

Assembly Insert new blister pack for next wear period.

- 1 Place the lower part of the clear plastic Blister Pack into the black holder.

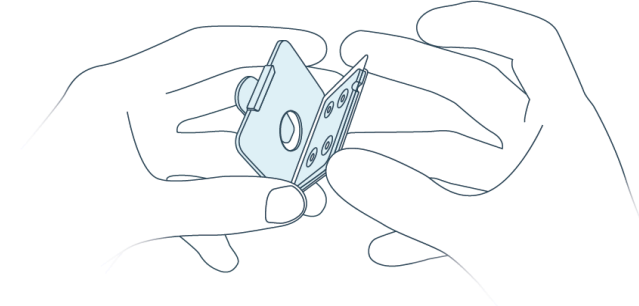


- 2 Press the top of the Blister Pack towards the holder until it snaps into place.

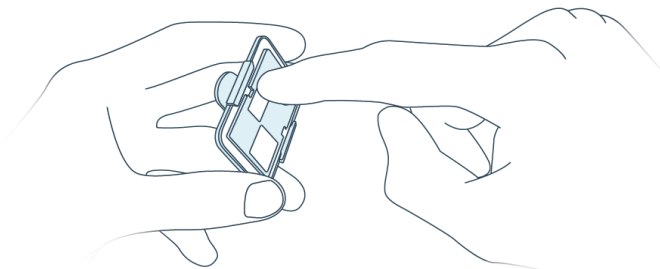


***WARNING:** Piercing, opening, tearing, or cutting the blister pack may cause permanent damage to the internal components of your dosimeter—rendering it unable to accurately and effectively measure dose. Damaged Blister Packs may incur a damage fee.

Disassembly Remove old Blister Pack (end of wear period) for processing.

- 1 Remove the Blister Pack by turning the clip (on the back of the holder) to expose the hole then pressing through with your finger.

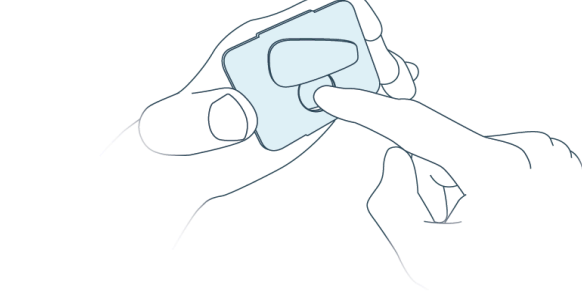

- 2 **Keep the black holder** and return the plastic Blister Pack for processing.

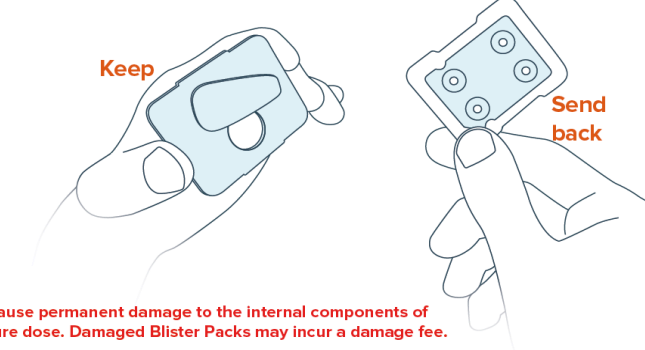

Problems with
a dosimeter
or shipment?

Contact Us

Customer Service

pds@radpro.co.nz

0508-RADPRO

radpro
Radiation Protection Services Limited